

Quick Start Guide

Product Codes: PB501, PC502, PV503, PU504



This guide provides a summary of the procedures for using Photopette® device from Tip Biosystems Pte Ltd. It is meant to be used only after a User has been fully trained according to the device's User Manual and lab procedures. This guide assumes that the User is fully aware of the health and safety issues associated with this system and its components as described in the system's User Manual which can be found on our company's website stated below.

Legal

Photopette® and CuveTip™ are trademarks of Tip Biosystems Pte Ltd

All goods and services are sold subject to the terms and conditions of sale of Tip Biosystems Sales Agreement. A copy of these terms and condition are available upon request.

©2016 Tip Biosystems Pte Ltd - All rights reserved.

First published March 2017

<http://www.tipbiosystems.com>

Version 1.0.2

1. Download the Photopette® App




2. Connecting to your Photopette®

- Switch on your Photopette® by pressing the button located on the side of your device.
- In the Photopette® App, navigate to “Devices”.
- Select your Photopette® from the list.

3. Selecting Application Types

- Select “Measurement Type” on the home screen.
- Select the intended application from the list. Note that applications are device specific.
- After selecting the desired application, press “Start Measurement”.
- **At this stage, ensure a CuveTip™ has been placed on your Photopette®.**

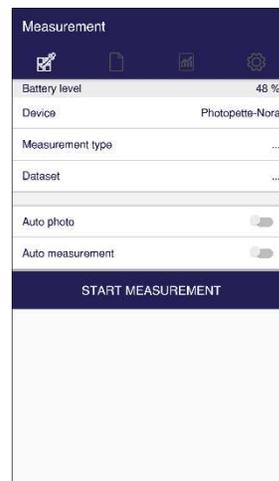

4. Measuring your reference sample

- Insert CuveTip™ into your reference sample (blank).
- The optimal depth of insertion is highlighted in the Figure on the right.
- Press the measurement button located on top of your Photopette®.



5. Measuring your sample

- Place a new CuveTip™ on your Photopette®.
- Insert CuveTip™ into your measurement sample.
- Press the measurement button.
- **Do not move the device when a measurement is in progress.**

6. Results

- Results will be displayed on the measurement screen.
- To review the data at a later point, navigate to “Dataset”.
- Here, you may also choose to export your data via Email or Google Drive.

λ (nm)	A	T (%)
600	2.528	0.296
600	2.694	0.202
600	2.360	0.436
600	2.276	0.530

Buttons: Press to initiate Measurement MEASUREMENT

Time	Measurement Type
02 Aug 17, 05:48:40 PM	Turbidity
03 Aug 17, 11:29:49 AM	ssRNA
03 Aug 17, 11:24:25 AM	ssDNA
03 Aug 17, 11:21:41 AM	ssDNA
03 Aug 17, 11:17:33 AM	ssRNA
03 Aug 17, 11:14:32 AM	ssRNA
03 Aug 17, 11:06:22 AM	ssDNA
03 Aug 17, 10:49:51 AM	

